



**Eldex**<sup>®</sup>

Eldex Laboratories, Inc.  
30 Executive Ct.  
Napa, CA 94558-6278

Tel: (707)224-8800  
(800)969-3533  
Fax: (707)224-0688

## Procedures

### Zero Micrometer

- 1) Loosen micrometer lock ring.
- 2) Turn on pump.
- 3) Back off micrometer to +.250.
- 4) Adjust micrometer inward using minimal force. Micrometer should move only when rocker arm is not in contact with micrometer face.
- 5) Continue this process until rocker arm is always against the micrometer face and no further adjustment is possible when only light pressure is used. This is the zero point.
- 6) Loosen the Allen set screw on the end of the micrometer.
- 7) Adjust micrometer so that it reads .000.
- 8) Tighten Allen set screw.